

# CHIP-IT® PBMC

(version A2)

Catalog No. 53042

## **Active Motif North America**

1914 Palomar Oaks Way, Suite 150  
Carlsbad, California 92008, USA  
Toll free: 877 222 9543  
Telephone: 760 431 1263  
Fax: 760 431 1351

## **Active Motif Europe**

Avenue Reine Astrid, 92  
B-1310 La Hulpe, Belgium  
UK Free Phone: 0800 169 31 47  
France Free Phone: 0800 90 99 79  
Germany Free Phone: 0800 181 99 10  
Telephone: +32 (0)2 653 0001  
Fax: +32 (0)2 653 0050

## **Active Motif Japan**

Azuma Bldg, 7th Floor  
2-21 Ageba-Cho, Shinjuku-Ku  
Tokyo, 162-0824, Japan  
Telephone: +81 3 5225 3638  
Fax: +81 3 5261 8733

Copyright 2014 Active Motif, Inc.

Information in this manual is subject to change without notice and does not constitute a commitment on the part of Active Motif, Inc. It is supplied on an “as is” basis without any warranty of any kind, either explicit or implied. Information may be changed or updated in this manual at any time.

This documentation may not be copied, transferred, reproduced, disclosed, or duplicated, in whole or in part, without the prior written consent of Active Motif, Inc. This documentation is proprietary information and protected by the copyright laws of the United States and international treaties.

The manufacturer of this documentation is Active Motif, Inc.

© 2014 Active Motif, Inc., 1914 Palomar Oaks Way, Suite 150; Carlsbad, CA 92008. All rights reserved.

All trademarks, trade names, service marks or logos referenced herein belong to their respective companies.

<b>TABLE OF CONTENTS</b>	<b>Page</b>
<b>Overview</b> .....	1
<b>Flow Chart of Process</b> .....	2
<b>Kit Performance and Benefits</b> .....	3
Protocol Overview and Time Table .....	4
<b>Kit Components and Storage</b>	
ChIP-IT PBMC Kit .....	5
Additional Materials Required .....	6
<b>Protocols – Experimental Set Up</b>	
Cell Recommendations .....	7
Buffer Preparation .....	7
Quick Chart for Preparing Buffers .....	8
Recommendations .....	9
<b>Protocols – Preparation of Sheared Chromatin</b>	
Section A. Cell Fixation .....	10
Section B. Chromatin Sonication .....	11
<b>Protocols – Chromatin Immunoprecipitation</b>	
Section C. Immunoprecipitation .....	14
<b>Protocols – ChIP DNA Purification</b>	
Section D. Reversal of Cross-links and DNA Purification .....	16
<b>Protocols – ChIP DNA Analysis</b>	
Section E. Quantitative PCR (qPCR) .....	17
Section F. ChIP-Seq .....	18
<b>Appendix</b>	
Section G. qPCR Primer Design and Data Analysis .....	20
Section H. Troubleshooting Guide .....	22
Section I. Related Products .....	24
<b>Technical Services</b> .....	26

## Overview

---

Chromatin Immunoprecipitation (ChIP) is a powerful tool for studying protein/DNA interactions, including transcription factors, co-regulatory proteins, modified histones, chromatin-modifying enzymes and polymerases because it enables identification of the localization of proteins bound to specific DNA loci. When used in combination with whole-genome analysis such as ChIP-Seq or ChIP-chip, insights are possible into gene regulation, gene expression, mechanisms of chromatin modification and pathway analysis.


However, ChIP can be technically demanding. The method requires high-quality antibodies to recognize the fixed, target-bound proteins of interest, and an efficient means to precipitate the antibody/chromatin complex (usually protein A or G beads). In addition, specialized buffers, inhibitor cocktails and blocking reagents are required to minimize non-specific enrichment and reduce protein degradation. The difficulties of ChIP are further compounded when working with difficult to lyse cells, such as peripheral blood mononuclear cells (PBMCs), including lymphocytes (T cells, B cells and NK cells) and monocytes.

Active Motif's ChIP-IT® PBMC Kit is designed to overcome these obstacles and generate high quality chromatin from PBMCs (including lymphocytes and monocytes) for analysis using the highly sensitive, low background ChIP reagents provided in the Kit. These optimized reagents and protocols translate into better ChIP data.

The ChIP-IT PBMC Kit contains sufficient reagents to perform 16 chromatin preparation and immunoprecipitation reactions. We recommend the use of Active Motif's ChIP-IT® qPCR Analysis Kit in combination with the ChIP-IT PBMC Kit for complete data analysis. To learn about available ChIP-IT® Control Kits, control qPCR primer sets, ChIP-Seq validated antibodies, or Active Motif's EpiShear™ sonication devices, please visit our website at [www.activemotif.com/chip](http://www.activemotif.com/chip).

<b>product</b>	<b>format</b>	<b>catalog no.</b>
ChIP-IT® PBMC	16 rxns	53042
ChIP-IT® qPCR Analysis Kit	10 rxns	53029

## Flow Chart of Process

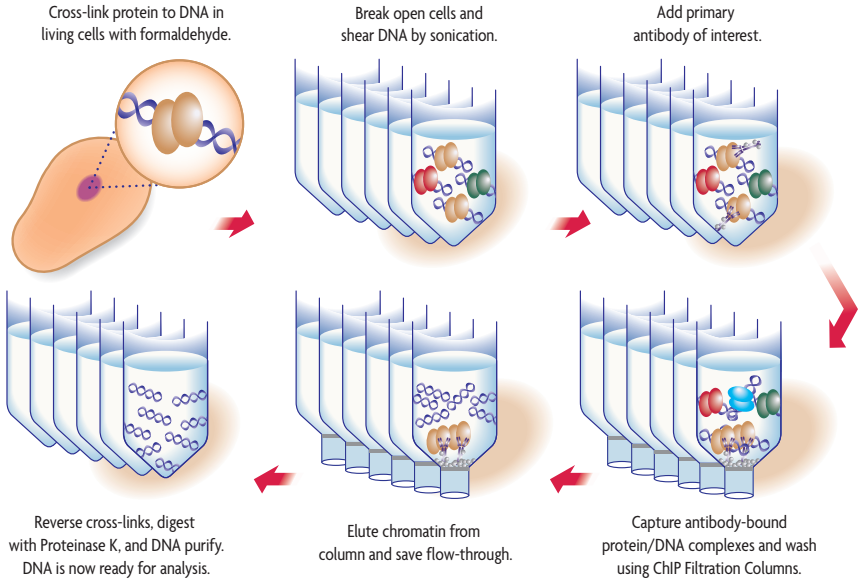


### Flow Chart of the ChIP-IT PBMC Assay.

In ChIP-IT PBMC, intact cells are fixed with formaldehyde, which cross-links and preserves protein/DNA interactions. DNA is then sheared into small fragments using sonication and incubated with an antibody directed against the DNA-binding protein of interest. The antibody-bound protein/DNA complexes are immunoprecipitated through the use of Protein G agarose beads and washed via gravity filtration. Following immunoprecipitation, cross-links are reversed, the proteins are removed by Proteinase K, and the DNA is recovered and purified. ChIP enriched DNA can be used for either gene-specific or whole-genome analysis.

## Kit Performance and Benefits


### ChIP-IT PBMC Advantages:

- Optimized buffers improve cell lysis and chromatin yields when working with PBMCs (including T cells, B cells, NK cells and monocytes)
- Highly sensitive DNA enrichment due to the inclusion of an antibody blocker which reduces background levels caused by non-specific binding events
- Filtration based washes are the easiest wash method available and result in increased consistency in multi-sample experiments
- Highly robust procedure has been validated in both qPCR and ChIP-Seq analysis

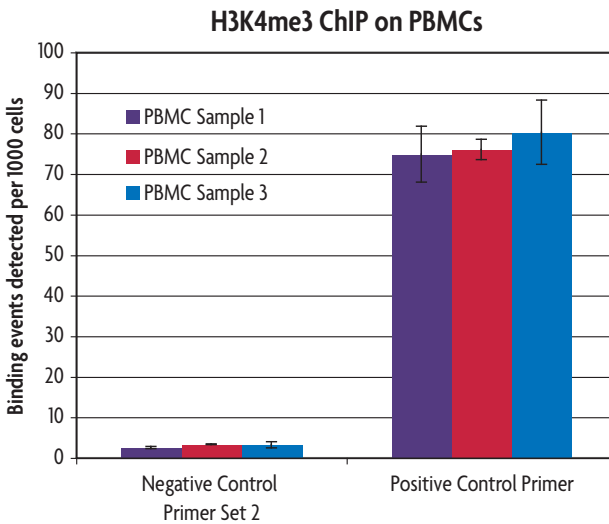

**Figure 1: Immunoprecipitation sensitivity of ChIP-IT High Sensitivity Kit.**

PBMC chromatin was prepared according to the manual using 20 million PBMC cells per reaction. ChIP was performed using the Histone H3K4me3 antibody on 10  $\mu$ g of chromatin. Following enrichment, qPCR was performed using the ChIP-IT qPCR Analysis Kit (Catalog No. 53029) to normalize the data. The negative control primer set (Catalog No. 71002) showed no enrichment, while the positive control primer sets showed nice enrichment of H3K4m3 in the three PBMC samples tested. The data represents triplicate values expressed as Binding events detected per 1,000 cells. To convert this scale to the percent of ChIP input recovered, divide the values by 1,000.

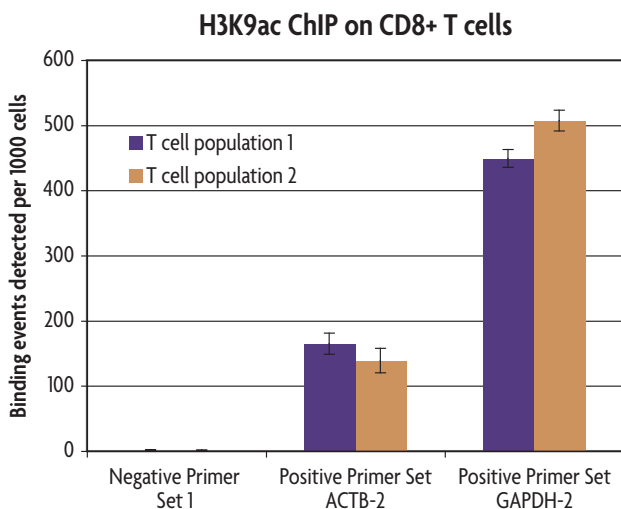

**Figure 2: ChIP-IT PBMC on CD8+ selected T cells.**

T cell chromatin was prepared according to the manual using 20 million CD8+ sorted T cells per reaction. ChIP was performed using the Histone H3K9ac antibody on 10 µg of chromatin. Following enrichment, qPCR was performed using the ChIP-IT qPCR Analysis Kit (Catalog No. 53029) to normalize the data. The negative control primer set (Catalog No. 71001) showed no enrichment, while both the ACTB-2 (Catalog No. 71005) and GAPDH-2 (Catalog No. 71006) positive control primer sets showed nice enrichment of H3K9ac in the two T cell populations tested. The data represents triplicate values expressed as Binding events detected per 1,000 cells. To convert this scale to the percent of ChIP input recovered, divide the values by 1,000.

## Protocol Overview and Time Table

	Required Time
Cell Fixation and Lysis	2 hours
Chromatin Sonication	40 minutes per sample
Assessment of Chromatin Size	Overnight incubation
Immunoprecipitation	Overnight incubation
Binding to Protein G agarose Beads	3 hours
Wash Immune Complexes	20 minutes
Reversal of Cross-links	2.5 hours
DNA Purification	15 minutes
qPCR Analysis	2 hours

## Kit Components and Storage

---

Please store each component at the temperature indicated in the table below. **Do not re-freeze the Protein G Agarose Beads. Once thawed, Protein G beads should be stored at 4°C.**

Reagents	Quantity	Storage
RNase A (10 µg/µl)	40 µl	-20°C
Proteinase K (10 µg/µl)	200 µl	-20°C
Blocker	100 µl	-20°C
100 mM PMSF	500 µl	-20°C
Protease Inhibitor Cocktail (PIC)	500 µl	-20°C
Precipitation Buffer	1.5 ml	-20°C
Carrier	35 µl	-20°C
10X PBS	120 ml	-20°C
Protein G Agarose Beads*	500 µl	4°C
Swelling Buffer	100 ml	4°C
TE pH 8.0	3 x 1.5 ml	RT
Detergent	25 ml	RT
Stop Solution	20 ml	RT
5 M NaCl	400 µl	RT
ChIP Filtration Columns	16 ea	RT
ChIP Buffer	35 ml	RT
Wash Buffer AM1	100 ml	RT
Elution Buffer AM4	2 x 1.5 ml	RT
DNA Purification Binding Buffer	50 ml	RT
3 M Sodium Acetate	500 µl	RT
DNA Purification Wash Buffer**	10 ml	RT
DNA Purification Elution Buffer	5 ml	RT
DNA Purification Columns	16 ea	RT

\* The Protein G Agarose Beads are shipped on dry ice and can be stored frozen until their first use. Once thawed, the Protein G beads **should not be re-frozen** by the customer. Protein G Agarose Beads should be stored at 4°C.

\*\*Requires the addition of ethanol before use.



## Additional materials required

- A ChIP-validated antibody directed against the protein of interest
- 37% formaldehyde solution with 10-15% methyl alcohol to prevent polymerization. Do not use paraformaldehyde
- 100% ethanol (absolute)
- 70% ethanol
- DNase-free H<sub>2</sub>O
- Centrifuge
- 15 and 50 ml conical tubes
- Metal spatula
- Dry ice for flash freezing
- Apparatus to rotate tubes end-to-end at 4°C (*e.g.* a Labquake from Barnstead/ThermoLyne with a tube holder for 1.5 ml microcentrifuge tubes or Glas-Col, Catalog No. 099A RD4512)
- Microcentrifuge (table top centrifuge 4°C) and microcentrifuge tubes
- 250 µl PCR tubes
- Thermocycler
- Phenol
- Chloroform/isoamyl alcohol (24:1)
- Spectrophotometer for DNA quantitation
- Pipettors and tips (filter tips are recommended)
- Sonicator (*e.g.* Active Motif's EpiShear™ Sonicator with a 1/8" probe (Catalog No. 53051) with the EpiShear™ Cooled Sonication Platform (Catalog No. 53080))
- Agarose gel electrophoresis apparatus
- (Optional) ChIP-IT® qPCR Analysis Kit (Catalog No. 53029)
- (Optional) Gene-specific qPCR primer pairs for enrichment analysis; see Appendix Section I
- (Optional) SYBR Green qPCR master mix (Bio-Rad Catalog No. 170-8882)

## Protocols – Experimental Set Up

---

**PLEASE READ THE ENTIRE PROTOCOL BEFORE STARTING!**

### Cell Recommendations

When planning an experiment, calculate the number of chromatin preparations you will require and determine the number of ChIP reactions you plan to perform on each chromatin preparation. Be sure to include the appropriate positive and negative control ChIP reactions in your calculations. Also, note that if you wish to analyze the effect of particular compounds on protein/DNA interactions, you should prepare chromatin from control (untreated) cells as a reference sample. PBMCs (including lymphocytes and monocytes) are difficult to lyse cells and as a result will yield approximately 30-50% as compared to cell lines. The minimum recommended number of cells that should be used for the preparation of chromatin is 10 million cells.

### Buffer Preparation

A Quick Chart for preparing the buffers needed for Section A: Cell Fixation is provided on page 8. Please refer to the descriptions below for complete details on buffer preparations.

#### Fixation Solution

Fixation solution should be prepared fresh before each experiment. For each cell pellet, prepare 5 ml of Fixation Solution by adding 0.5 ml 10X PBS to 4.36 ml sterile water in a 15 ml conical tube. Using appropriate precautions (*i.e.* safety glasses, gloves and lab coat), add 140  $\mu$ l 37% formaldehyde to the tube and vortex to mix.

#### Stop Solution

Is provided ready to use. Use 250  $\mu$ l per 5 ml Fixation Solution.

#### 100 mM PMSF and Protease Inhibitor Cocktail (PIC)

Thaw the PMSF and the PIC at room temperature until fully dissolved (approximately 30 minutes). Vortex gently and spin down briefly before use, then add to the buffers immediately before use.

#### Detergent

Is supplied ready to use.

#### PBS Wash Buffer

Prepare 10 ml PBS Wash Buffer for every sample. To a 15 ml conical tube add 8.5 ml sterile water, 1 ml 10X PBS, 0.5 ml Detergent and 20  $\mu$ l 100 mM PMSF. Mix by inverting. Place PBS Wash Buffer on ice to chill. PBS Wash Buffer can be prepared in large quantities and stored at 4°C for 6 months.

#### Swelling Buffer

Is supplied ready to use.

#### ChIP Buffer

Is supplied ready to use.

### DNA Purification Wash Buffer

The DNA Purification Wash Buffer requires the addition of ethanol before use. The final concentration of ethanol should be 80%. Add 40 ml of fresh 100% ethanol to the DNA Purification Wash Buffer bottle. Invert repeatedly. The buffer can be stored at room temperature after the addition of ethanol. The ethanol only needs to be added before the first use, after that the Wash Buffer is ready for use.

### 3M Sodium Acetate

It is important to check the sodium acetate before use to ensure that the salts have not precipitated out of solution. Once the sodium acetate is in solution it should be stored at room temperature.

### Protein G Agarose Beads

The supplied agarose beads require washing prior to use. Follow the instructions in the manual to wash the beads for use in the ChIP reactions. There is no need to pre-block the beads or pre-clear the sample. For best results, gently shake and invert the tube to resuspend the agarose beads. The beads settle quickly, and therefore should be resuspended just before pipetting. We recommend cutting 2 mm from the end of a pipet tip prior to pipetting to prevent the tip from becoming clogged. **Protein G Agarose Beads are shipped on dry ice and can be stored frozen until their first use. Once thawed, beads should not be re-frozen by the customer. Protein G Agarose Beads should be stored at 4°C.**

## Quick Chart for Preparing Buffers for Section A: Cell Fixation

Reagents to prepare	Components	Per cell pellet	4 cell pellets
Fixation Solution	Sterile H <sub>2</sub> O	4.36 ml	17.44 ml
	10X PBS	0.5 ml	2 ml
	37% Formaldehyde	140 µl	560 µl
	<b>TOTAL REQUIRED</b>	<b>5 ml</b>	<b>20 ml</b>
Stop Solution	<b>TOTAL REQUIRED</b>	<b>250 µl</b>	<b>1 ml</b>
PBS Wash Buffer	Sterile H <sub>2</sub> O	8.5 ml	34 ml
	10X PBS	1 ml	4 ml
	Detergent	0.5 ml	2 ml
	100 mM PMSF	20 µl	80 µl
	<b>TOTAL REQUIRED</b>	<b>10 ml</b>	<b>40 ml</b>
Swelling Buffer	Swelling Buffer	5 ml	20 ml
	Protease Inhibitor Cocktail	5 µl	20 µl
	100 mM PMSF	5 µl	20 µl
	<b>TOTAL REQUIRED</b>	<b>5 ml</b>	<b>20 ml</b>
Detergent	<b>TOTAL REQUIRED</b>	<b>375 µl</b>	<b>1.5 ml</b>
ChIP Buffer	ChIP Buffer	500 µl	2 ml
	Protease Inhibitor Cocktail	5 µl	20 µl
	100 mM PMSF	5 µl	20 µl
	<b>TOTAL REQUIRED</b>	<b>510 µl</b>	<b>2 ml</b>

## Recommendations

### ChIP-validated Antibody

We recommend using 4 µg antibody per ChIP reaction in a maximum volume of 30 µl. However, this will vary according to the affinity of the antibody and the quality of the chromatin; you may need to use more of a particular antibody. ChIP antibodies must recognize fixed, native protein that is bound to DNA and/or complexed with other proteins. Many antibodies that perform well in other applications do not perform in ChIP. Thus, ChIP performed with an antibody that has not been ChIP-validated must include appropriate controls (such as Active Motif's ChIP-IT Control qPCR Kits, Catalog Nos. 53026, 53027 and 53028) to validate the chromatin preparation and the ChIP methodology. To see a list of available ChIP-validated antibodies available from Active Motif, please visit [www.activemotif.com/chipabs](http://www.activemotif.com/chipabs).

### Chromatin Shearing Tips

We suggest using a probe sonicator (*i.e.* Active Motif's EpiShear Probe Sonicator) which employs a direct sonication method to prepare chromatin for use in the ChIP-IT PBMC Kit. Indirect sonication systems may require longer sonication times to achieve optimal chromatin shearing. ChIP experiments usually require chromatin that has been sheared to a size of 200-1200 bp. In general, shearing efficiency is improved through the use of a small shearing volume and a V-bottom tube rather than a round-bottom tube. Also, note that shearing is inefficient if the chromatin sample becomes emulsified with air bubbles. To determine the appropriate shearing level for your sample, set up a "practice" tube containing only ChIP Buffer. Slowly increase the sonication amplitude until foaming starts to occur. Reduce the amplitude setting down slightly and mark this as the highest possible intensity to use without foaming. If a chromatin preparation becomes emulsified inadvertently, discontinue shearing and centrifuge the sample at maximum speed for 4 minutes at 4°C in a microcentrifuge to remove trapped air. Finally, to prevent overheating and denaturation of chromatin, samples should be kept on ice as much as possible during shearing, and shearing should be performed discontinuously (*i.e.* sonicate for 20 seconds, then place on ice/water for 30 seconds, sonicate again for 20 seconds, *etc.*). If possible, shear while on ice or use Active Motif's EpiShear™ Cooled Sonication Platform (Catalog No. 53080) to help regulate sample temperature.

### Chromatin Quantity

It is recommended to use 10-30 µg chromatin per IP reaction. However, if chromatin is in limited supply lower amounts may be used due to the high quality of the DNA obtained from the ChIP-IT PBMC Kit. If ChIP-Seq will be performed following enrichment, please read the recommendations listed in Section F prior to starting the protocol.

### Safety Precautions

Formaldehyde and PMSF are highly toxic chemicals. Appropriate safety precautions (*i.e.* safety glasses, gloves and lab coat) should be used. Also, formaldehyde is highly toxic by inhalation and should be used only in a ventilated hood. Finally, chromatin sonication should be performed in a biosafety hood if the chromatin is extracted from biohazardous or infectious materials.

# Protocols – Preparation of Sheared Chromatin

---

## Section A: Cell Fixation

We recommend using a minimum of 10 million cells for the preparation of chromatin.

1. Collect your PBMCs, T cells, B cells, NK cells or monocytes according to your preferred protocol. Transfer the appropriate number of cells required for each chromatin preparation into a conical tube and centrifuge at  $1,250 \times g$  at  $4^{\circ}\text{C}$  to collect the cells. If there are cells that are floating or are stuck to the sides of the tube following the spin, use a metal spatula to scrape the cells down into solution and repeat the centrifugation at  $1,250 \times g$  at  $4^{\circ}\text{C}$  for 5 minutes. Continue to repeat the scraping and spin as many times as needed to pellet all the cells. Remove and discard supernatant.

**Note:** PBMCs may stick to certain plastics. Therefore, it is recommended to use a metal spatula for cell scraping to avoid losing any sample material.

2. Flash freeze the cell pellet(s) by immersing conical tube into dry ice. Incubate on dry ice for 10 minutes. Cell pellets may be stored at  $-80^{\circ}\text{C}$  at this stage, or proceed with the next step.
3. Prepare fresh Fixation Solution. To fix cells, add 5 ml of freshly prepared Fixative Solution to the cell pellet and resuspend cells by pipetting up and down. Incubate at room temperature for 15 minutes on a roller.
4. Stop the fixation reaction by adding 250  $\mu\text{l}$  Stop Solution to the existing fixation solution. Swirl to mix and incubate at room temperature for 5 minutes on a roller.
5. Following the incubation, add 250  $\mu\text{l}$  Detergent to each conical tube. Invert to mix. Pellet the cells by centrifugation for 5 minutes at  $1,250 \times g$  at  $4^{\circ}\text{C}$ .
6. If there are cells that are floating or are stuck to the sides of the tube following the spin, use a metal spatula to scrape the cells down into solution and repeat the centrifugation at  $1,250 \times g$  at  $4^{\circ}\text{C}$  for 5 minutes. Continue to repeat the scraping and spin as many times as needed to pellet all the cells.
7. Remove the supernatant and discard. Resuspend the pellet(s) in 5 ml ice-cold PBS Wash Buffer by pipetting up and down. Keep samples ice-cold for the remainder of the procedure.
8. Centrifuge for 5 minutes at  $3,200 \times g$  at  $4^{\circ}\text{C}$ . The cell pellet will tend to be loose, so carefully remove the supernatant and discard taking care to avoid disturbing the pellet. Wash the pellet(s) a second time in 5 ml ice-cold PBS Wash Buffer by pipetting up and down. Centrifuge for 5 minutes at  $3,200 \times g$  at  $4^{\circ}\text{C}$ . Carefully remove the supernatant and discard.
9. Flash freeze the cell pellet(s) by immersing conical tube into dry ice. The freeze/thaw will help facilitate cell lysis. Incubate on dry ice for 10 minutes. Cell pellets may be stored at  $-80^{\circ}\text{C}$  at this stage, or proceed with the next step.
10. Resuspend each pellet(s) in 5 ml ice-cold Swelling Buffer supplemented with 5  $\mu\text{l}$  PIC and 5  $\mu\text{l}$  100 mM PMSF. Pipet up and down to mix.
11. Incubate on ice for 30 minutes.

12. Add 125 µl Detergent to each cell pellet. Vortex for 30 sec on highest setting to lyse the cells.

**Note:** We do not recommend the use of a dounce homogenizer to assist in cell lysis as the transfer process to and from the dounce increases sample loss. If cells continue to have difficulty lysing, we suggest additional sonication.

13. Centrifuge for 10 minutes at 3,200 x g at 4°C. Remove and discard the supernatant.
14. Resuspend each pellet in 500 µl ChIP Buffer supplemented with 5 µl PIC and 5 µl 100 mM PMSF. Transfer the contents to a new 2 ml microcentrifuge tube.
15. Incubate on ice for 10 minutes. Proceed to Step B: Chromatin Sonication.

## Section B. Chromatin Sonication

The section below describes the fragmentation of chromatin using sonication. Sonication results may vary depending on cell type and sonication device being used. This protocol has been validated using Active Motif's EpiShear™ Probe Sonicator in combination with an EpiShear™ Cooled Sonication Platform to maintain probe height and temperature consistency between samples. We do not recommend sonication of samples containing less than  $1 \times 10^7$  cells and/or 350 µl volume.

The ChIP Buffer has been optimized for immunoprecipitation performance, however, due to its unique composition optimization of sonication conditions may be required. To maintain the high sensitivity of the assay, we recommend using our buffer system and altering the sonication time and/or amplitude of your sonication system to achieve the desired fragmentation (*e.g.* some systems may require as much as a three-fold increase in sonication time to improve chromatin shearing). Please pay particular attention to our protocol regarding the processing of input chromatin for agarose gel analysis prior to the chromatin immunoprecipitation reaction as many steps may differ from traditional ChIP protocols and failure to follow the outlined procedure may lead to artifacts in the gel images as shown in Figure 3 on page 13.

1. Place the 2 ml microcentrifuge tube containing the chromatin into the tube cooler or packed ice. Open cap and submerge the microtip into the liquid until the microtip is approximately 5 mm from the bottom of the tube. Sonicate according to optimized settings for the cell type being used (see Recommendations on page 9). A recommended starting range for PBMCs is: 4 rounds of sonication at 42% amplitude, pulse for 30 seconds on and 30 seconds off for a total sonication “on” time of 5 minutes per round. (Total sonication “on” time is 20 minutes and total elapsed time is 40 minutes per sample.)
2. Spin tubes at 4°C in a microcentrifuge at maximum speed for 5 minutes to pellet the cellular debris.
3. Transfer 25 µl of each chromatin preparation into a 250 µl PCR tube for analysis of shearing efficiency and chromatin quantification. This sample will be used to generate the Input DNA.
4. Aliquot the remainder of each chromatin preparation into 1.5 ml microcentrifuge tubes. We recommend making aliquots of 250 µl volume.
5. Flash freeze the chromatin by immersing tubes into dry ice for 10 minutes and store at -80°C.


**Note:** The size of the chromatin sonication should be verified before proceeding to the immunoprecipitation step.

### *Input Preparation*

6. To each 25  $\mu$ l chromatin preparation from Step 5 above, add 175  $\mu$ l TE pH 8.0 and 2  $\mu$ l RNase A. Cap the PCR tubes and vortex to mix.
7. Incubate in a thermocycler at 37°C for 1 hour.
8. Add 5  $\mu$ l Proteinase K to each tube, vortex and incubate in a thermocycler at 37°C for 3 hours.
9. Add 10  $\mu$ l 5 M NaCl, vortex and incubate at 65°C for 6-16 hours to reverse cross-links.
10. Remove tubes from the thermocycler and add 250  $\mu$ l phenol and 125  $\mu$ l chloroform:isoamyl alcohol (24:1). Vortex vigorously and spin tubes in a room temperature microcentrifuge at maximum speed for 2 minutes.
11. Transfer each upper aqueous layer to a new 1.5 ml microcentrifuge tube and add 250  $\mu$ l chloroform:isoamyl alcohol (24:1). Vortex vigorously and spin tubes in a room temperature microcentrifuge at maximum speed for 2 minutes.
12. Transfer the upper aqueous layer to a new 1.5 ml microcentrifuge tube. Add 83  $\mu$ l Precipitation Buffer, 2  $\mu$ l Carrier and 900  $\mu$ l absolute ethanol. Vortex to mix and chill at -80°C for 30 minutes to overnight.
13. Spin at 4°C in a microcentrifuge at maximum speed for 15 minutes.
14. Carefully remove the supernatant taking care not to disturb the pellet. Wash the pellet with 500  $\mu$ l 70% ethanol and spin at 4°C in a microcentrifuge at maximum speed for 5 minutes.
15. Carefully remove the supernatant taking care not to disturb the pellet. Remove residual ethanol with a pipet tip. Leave the tubes uncapped and air dry for 10-15 minutes.
16. When the pellets are dry, add 25  $\mu$ l DNA Purification Elution Buffer to each tube. Incubate at room temperature for 10 minutes. Then vortex to ensure the pellet is completely resuspended. This solution contains your Input DNA.
17. Read the absorbance of each sample on a NanoDrop or other spectrophotometer at 260 nm to determine the DNA concentration of each chromatin preparation. Set aside 500 ng of DNA for analysis as described in Step 18. Store the remaining Input DNA at -20°C.
18. Analyze each chromatin preparation on an agarose gel by following the instructions below.
  - a. Prepare 500 mM NaCl by adding 2  $\mu$ l 5M NaCl to 18  $\mu$ l sterile water. Vortex to mix.
  - b. Transfer 500 ng of Input DNA to a 250  $\mu$ l PCR tube and add 1  $\mu$ l 500 mM NaCl. Adjust the final volume to between 10  $\mu$ l with sterile water if needed.
  - c. Heat samples in a thermocycler at 100°C for 20 minutes followed by ramping the temperature down to 50°C.
  - d. Remove tubes from the thermocycler and incubate at room temperature for 5 minutes.
  - e. Add gel loading buffer to each sample and run on a 1.5% agarose gel. Include 100 bp

and 1 kb DNA ladders to analyze chromatin size. DNA should appear as a smear anywhere between 200-1200 bp.

19. If chromatin preparations were successful, the aliquots stored at  $-80^{\circ}\text{C}$  from Section B, Step 5 can be used to perform the ChIP reactions in Section C.

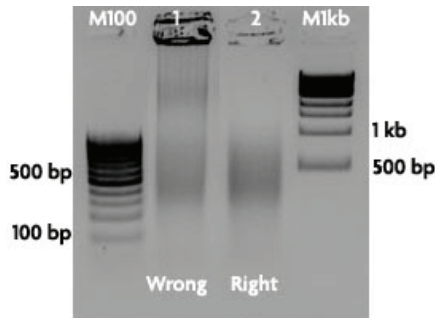

**Figure 3: Validation of chromatin shearing efficiency following reversal of cross-links overnight at  $65^{\circ}\text{C}$ .**

Chromatin preparations of MCF-7 cells were fixed and sonicated using the EpiShear™ Probe Sonicator and EpiShear™ Cooled Sonication Platform from Active Motif. Input DNA was prepared in duplicate according to Section B, Steps 6-17 in the manual. In Sample 1, Step 18 was not performed and 500 ng of input DNA was loaded directly onto a 1.5% agarose gel without receiving the addition of NaCl and incubation at  $100^{\circ}\text{C}$ . The omission of Step 18 has caused a buffer artifact that makes the DNA appear larger on a gel. The duplicate sample, Sample 2, was processed according to the manual instructions and included the addition of NaCl and incubation at  $100^{\circ}\text{C}$  as stated in Step 18. Analysis of 500 ng of this input DNA on a 1.5% agarose gel shows the expected fragmentation between 200-1200 bp. The difference in DNA sizing on the gel between the two samples illustrates the importance of following the protocol recommendations regarding the processing of input chromatin for agarose gel analysis prior to chromatin immunoprecipitation. Omission of key steps can lead to inaccurate analysis of chromatin shearing efficiency. If the protocol steps were followed and the DNA fragments fall outside of the recommended range, sonication conditions should be further optimized.



# Protocols – Chromatin Immunoprecipitation

## Section C. Immunoprecipitation

Successful chromatin immunoprecipitation depends on the quality of the ChIP antibody and the abundance of the target protein. We recommend using 10 – 30 µg chromatin per IP reaction, although less can be used if chromatin yields are limiting.

1. Thaw sonicated chromatin on ice. Spin chromatin at 4°C in a microcentrifuge at maximum speed for 2 minutes.
2. Set up the ChIP reactions by adding the components in the order shown in Table 1 below to 1.5 ml microcentrifuge tubes. Be sure to use the DNA concentration that was determined for your sonicated chromatin sample to calculate the volume to use.
3. In a 1.5 ml microcentrifuge tube prepare the antibodies to be used in the ChIP reactions. Use a separate tube for each antibody. To the tube add 5 µl Blocker and 4 µg ChIP antibody. (Antibody volume should not exceed 30 µl per reaction). Incubate Antibody/Blocker mix for 1 minute at room temperature and then add to the ChIP reactions.

**Table 1**

Reagent	1 reaction
Sheared Chromatin (10 - 30 µg)	X µl
ChIP Buffer	adjust up to 200 µl
Protease Inhibitor Cocktail (PIC)	5 µl
Antibody/Blocker mix (from Step 3)	not to exceed 35 µl
<b>Maximum Volume Allowed</b>	<b>240 µl</b>

**Note:** If your chromatin is too dilute to use within a 200 µl volume, we suggest preparing two immunoprecipitation reactions each using 200 µl sheared chromatin and then pooling the IP reactions together when loading sample onto the ChIP filtration column in Section C, Step 11.


4. Cap tubes and incubate on an end-to-end rotator overnight at 4°C.
5. The Protein G agarose beads require washing before use. Transfer 30 µl Protein G agarose beads for each IP reaction to a 1.5 ml microcentrifuge tube. Add an equal volume of TE, pH 8.0 and invert to mix. Spin at 1250 x g in a microcentrifuge for 1 minute. Remove the supernatant equivalent to the volume of TE added to the agarose beads.

**Note:** Before pipetting the Protein G agarose beads, they should be fully resuspended by inverting the tube. When pipetting the beads, cut 2 mm from the end of a pipet tip to prevent the tip from becoming clogged.

6. Wash the beads a second time with the same volume of TE, pH 8.0. Invert to mix. Spin at 1250 x g for 1 minute in a microcentrifuge. Remove the supernatant equivalent to the volume

of TE added to the agarose beads. The beads are now ready to use.

7. Spin the ChIP reactions at 1250 x g for 1 minute to collect liquid from the inside of the cap.
8. Using a cut pipet tip, add 30  $\mu$ l washed Protein G agarose beads to each immunoprecipitation reaction. Cap tubes and incubate on an end-to-end rotator at 4°C for 3 hours.
9. Label a ChIP Filtration Column for each ChIP reaction. Remove the tab from the bottom of the column and place in an empty 1 ml pipet tip box as a holder (see Figure 4 below).
10. Remove ChIP reactions from rotator and spin at 1250 x g for 1 minute to collect liquid from inside of the cap.
11. Add 600  $\mu$ l ChIP Buffer to each ChIP reaction then transfer the entire reaction (including the protein G agarose beads) to its labeled column. Allow flow-through to occur by gravity.
12. During the gravity flow, transfer 100  $\mu$ l per ChIP reaction of Elution Buffer AM4 to a 1.5 ml microcentrifuge tube and allow to pre-warm at 37°C during the wash steps.
13. Wash each column with 900  $\mu$ l Wash Buffer AM1. Let stand for 3 min.
14. Repeat Step 13 four more times for a total of five washes.
15. Transfer columns to a new 1.5 ml microcentrifuge tube and spin in a room temperature microcentrifuge at 1250 x g for 3 minutes to remove residual Wash Buffer.
16. Following the spin, transfer the ChIP Filtration Columns to new 1.5 ml microcentrifuge tubes. Add 50  $\mu$ l 37°C Elution Buffer AM4 to each column. Incubate at room temperature for 5 minutes. Spin in a room temperature microcentrifuge at 1250 x g for 3 minutes.
17. With columns remaining in the same microcentrifuge tube, add another 50  $\mu$ l 37°C Elution Buffer AM4 to each column. Incubate at room temperature for 5 minutes and spin in a room temperature microcentrifuge at 1250 x g for 3 minutes.
18. Discard the ChIP Filtration Columns. The flow-through (~100  $\mu$ l volume) contains the ChIP DNA. Proceed to Section D: DNA purification.

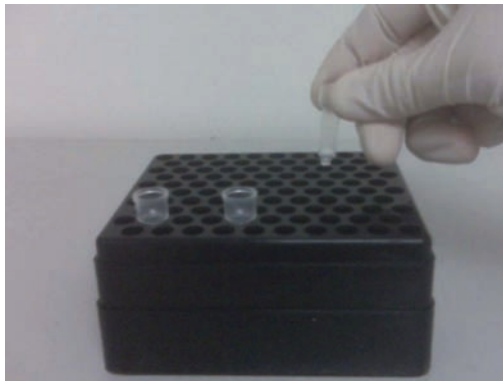

**Figure 4: Using the ChIP Filtration Columns.**

Remove the tab from the bottom of the ChIP Filtration Columns and place columns in an empty 1 ml pipet tip box to perform the wash steps.

### Section D. Reversal of Cross-links and DNA Purification

1. Transfer each eluted ChIP DNA to a 250  $\mu$ l PCR tube and add 2  $\mu$ l Proteinase K. Vortex to mix and heat in a thermocycler at 55°C for 30 minutes and then increase the temperature to 80°C for 2 hours.
2. Transfer the DNA to a 1.5 ml microcentrifuge tube and add 5 volumes (500  $\mu$ l) DNA Purification Binding Buffer to each tube and vortex to mix. Adjust the pH with 5  $\mu$ l 3M Sodium Acetate. The sample should be bright yellow in color to indicate a proper pH. If your sample is not bright yellow, please refer to the Troubleshooting guide in the Appendix for details to adjust pH prior to loading the sample into the purification column.
3. For each sample, place a DNA purification column (AM #103928) in the collection tube and add each pH adjusted sample to its own column. Close the cap on each column, place them with the collection tubes in a microcentrifuge and spin them at 14,000 rpm for 1 minute.
4. Remove the column from the collection tube, then remove and discard the flow through from the collection tube. Return the column to the collection tube.
5. Prepare DNA Purification Wash Buffer (AM #103497) before the first use. Follow the instruction on page 8 for the addition of ethanol prior to using the solution. Add 750  $\mu$ l DNA Purification Wash Buffer to each column and cap the column.
6. Spin at 14,000 rpm for 1 minute in a microcentrifuge.
7. Remove the column from the collection tube, then remove and discard the flow through from the collection tube. Return the column to the collection tube.
8. With the column cap open, spin at 14,000 rpm for 2 minutes in a microcentrifuge to remove any residual Wash Buffer from the column.
9. Transfer the column to a clean microcentrifuge tube. Depending on the method to be used for downstream analysis of the ChIP enriched DNA, use the appropriate elution volume. For ChIP-Seq or ChIP-chip applications, we recommend pre-warming the required elution buffer volume at 37°C for 5 minutes prior to use.
  - a. **For qPCR analysis:** add 100  $\mu$ l of DNA Purification Elution Buffer (AM #103498) to the center of the column matrix and incubate for 1 minute at room temperature. Spin at 14,000 rpm for 1 minute in a microcentrifuge. Add an additional 100  $\mu$ l DNA Purification Elution Buffer to the column and incubate for 1 minute at room temperature. Spin at 14,000 rpm for 1 minute in a microcentrifuge. Total elution volume is 200  $\mu$ l.
  - b. **For ChIP-Seq or ChIP-chip analysis:** add 36  $\mu$ l of 37°C DNA Purification Elution Buffer (AM #103498) to the center of the column matrix and incubate for 1 minute at room temperature. Spin at 14,000 rpm for 1 minute in a microcentrifuge.
10. Discard column. Purified DNA may be stored at -20°C for future use.

## Protocols – ChIP DNA Analysis

---

### Section E. Quantitative PCR (qPCR)

ChIP DNA can be analyzed on a gene-specific basis using quantitative PCR (qPCR). Positive control and negative control PCR primer pairs should be included in every analysis to determine the fold enrichment. Negative control primers will amplify a region of the genome not bound by the antibody target of interest. Active Motif recommends the use of its ChIP-IT® qPCR Analysis Kit (Catalog No. 53029) for the analysis of qPCR data. The ChIP-IT qPCR Analysis Kit contains positive and negative control primer pairs, standard curve DNA, standard curve primers and a qPCR Analysis spreadsheet to perform the analysis calculations. Active Motif's analysis strategy determines primer efficiencies and the ChIP sample values are normalized according to input, primer efficiency, chromatin amount used in the ChIP reaction and resuspension volume. The ChIP-IT qPCR Analysis Kit provides consistency in data analysis and allows direct comparison across samples and experiments. If not using the ChIP-IT qPCR Analysis Kit, qPCR data normalization and graphing can be done using the methods described in Section I.

1. Below is an example qPCR reaction. Please follow the specific instructions for your qPCR instrument. We recommend using a commercially available SYBR Green qPCR master mix (e.g. Bio-Rad Cat # 170-8882) and preparing triplicate reactions.

Reagent	20 µl PCR reactions
2X SYBR Green master mix	10 µl
Primer mix (2.5 µM each primer)*	4 µl
Sterile water	1 µl
DNA sample (ChIP or Input)	5 µl
<b>Total volume</b>	<b>20 µl</b>

\* We recommend designing primers to perform at an annealing temperature of 58°C so that all qPCR reactions can be performed under identical conditions. An amplicon length of 75-150 bp is recommended. Active Motif offers validated species-specific qPCR primers designed according to these recommendations at [www.activemotif.com/chipprimers](http://www.activemotif.com/chipprimers).

2. Place tubes in a real time PCR instrument and program as below:  
95°C for 2 minutes  
(95°C for 15 seconds, 58°C for 20 seconds, 72°C for 20 seconds) for 40 cycles
3. If using the ChIP-IT qPCR Analysis Kit, please refer to the product manual for analysis instructions. Otherwise, follow the recommendations in Section G.

**Notes:** The polymerase chain reaction (PCR) process for amplifying nucleic acid is covered by U.S. Patent Nos. 4,683,195 and 4,683,202 assigned to Hoffmann-La Roche. Patents pending in other countries.

## Section F. CHIP-Seq

The CHIP-IT PBMC Kit has been extensively validated for whole genome analysis using CHIP-Seq. This process involves the preparation of libraries from CHIP DNA by the addition of adapter sequences to the ends of the DNA fragments. The library is then PCR amplified and validated prior to sequencing.

Commercial kits are available to prepare the CHIP DNA libraries for sequencing such as Illumina® CHIP-Seq DNA Sample Prep Kit (Catalog No. IP-102-1001) which is compatible with sequencing on HiSeq 2000, HiSeq 1000, HiScanSQ or Genome Analyzer. For alternative sequencing platforms, please select an appropriate library preparation method compatible with the sequencing apparatus to be used.

### A. General Recommendations

- We recommend using 30 µg chromatin for the CHIP Reactions, although less may be used if chromatin amounts are limiting or you are enriching for a highly abundant histone protein
- Library generation usually requires 10 ng of CHIP enriched DNA. However, by using the CHIP-IT PBMC Kit the enriched DNA is of such a high quality that lower amounts can be used. (See B. CHIP-Seq Requirements below for more details)
- Sequence library yields should be in the range of 1-2 micrograms
- 30 million sequencing reads is sufficient for most transcription factor and histone modifications
- 50 bp single end reads are sufficient for unique mapping and good CHIP-Seq data, although longer reads can be used
- Input DNA should be sequenced as a control reaction in order to identify false “peaks” and also to reveal regions of the genome that have been duplicated. Subtracting the input peaks from the experimental peaks will help to eliminate false data. Use 50 ng Input DNA (from Section B Step 17) for library generation.

### B. CHIP-Seq Requirements


It is commonly recommended that library generation of CHIP DNA requires 10 ng of CHIP enriched DNA. Obtaining 10 ng of CHIP DNA should be attainable when using good antibodies against abundant histone modifications. However, 10 ng is an unrealistic number for a sequence specific DNA binding transcription factor. Using the CHIP-IT PBMC Kit, we have routinely generated high quality libraries from sub-nanogram levels of CHIP DNA due to the high quality of the DNA obtained. For good CHIP-Seq data, quality enrichments with low background is more important than the total quantity of DNA recovered. Therefore, we recommend performing qPCR on known binding sites to verify enrichment levels against a negative control primer set using Active Motif's CHIP-IT qPCR Analysis Kit (Catalog No. 53029) to confirm DNA quality rather than quantifying the DNA.

### C. qPCR prior to CHIP-Seq

To perform qPCR for verification of the quality of the enriched CHIP DNA, follow the instructions for qPCR in Section E with the following modification. The CHIP DNA will need to be diluted prior

to use in qPCR due to the reduced elution volume. Dilute 6 µl of ChIP DNA in 94 µl DNA Purification Elution Buffer (AM #103498). Use 5 µl of the diluted ChIP DNA per qPCR reaction.

Quality DNA that is suitable for use in ChIP-Seq should show enrichment of known binding sites over negative control primer sets of 5-fold (See Figure 5 below). If the qPCR enrichments are satisfactory then the remaining 30 µl of ChIP DNA can be used for library generation.

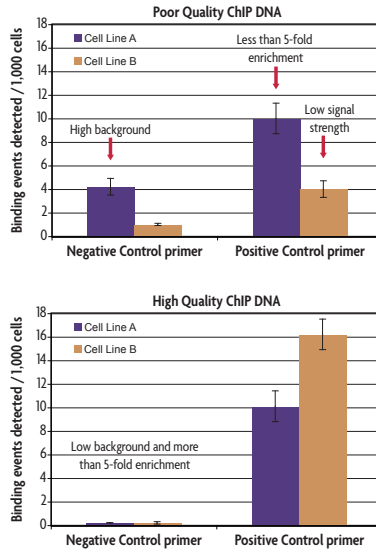

**Figure 5: Comparison of qPCR results showing good versus poor enrichment over negative control primers.**

Data shows qPCR results analyzed using the ChIP-IT qPCR Analysis Kit with the Human Negative Control Primer as a reference and a gene-specific positive control primer. In the top image the negative control primer set for Cell Line A gives high background levels with Binding events detected per 1,000 cells above a value of 2, while Cell Line B has positive control signal levels below 5 Binding events detected per 1,000 cells. The bottom image shows low background in the negative control primer set with Binding events detected per 1,000 cells below a value of 2. The fold enrichment of the positive control primer set exceeds 5-fold. Only the ChIP DNA from the bottom sample is recommended for use in ChIP-Seq.

#### D. Library Generation

ChIP DNA from the ChIP-IT PBMC Kit has been extensively validated for ChIP-Seq using the Illumina® platform. ChIP DNA can be used to generate single or paired-end libraries using the standard Illumina® library protocols.

To download an up to date version of the Illumina® library generation protocols, please visit the Illumina® website at [http://support.illumina.com/sequencing/sequencing\\_kits/chip-seq\\_dna\\_sample\\_prep\\_kit.ilmn](http://support.illumina.com/sequencing/sequencing_kits/chip-seq_dna_sample_prep_kit.ilmn), or search for Catalog No. IP-102-1001.

### Section G: qPCR Primer Design and Data Analysis

#### A. Design of the primers

- Design and analyze your potential primer pairs using an *in silico* PCR program (i.e. Primer3 at <http://frodo.wi.mit.edu/> or the UCSC Genome Browser at <http://genome.ucsc.edu/cgi-bin/hgPcr>).
- Primers that dimerize should be avoided, as they will be bound by SYBR Green, which will compromise accurate quantitation. You can test your primers for self-complementarity and secondary structure at [http://frodo.wi.mit.edu/cgi-bin/primer3/primer3\\_www.cgi](http://frodo.wi.mit.edu/cgi-bin/primer3/primer3_www.cgi).
- Ideally, the amplicons should be 75-150 bp in length.
- For use with the ChIP-IT qPCR Analysis Kit, primers should be designed to anneal optimally at 58°C with a recommended length of 18-22 bp.
- Active Motif offers ChIP Control qPCR primer sets validated to work in our ChIP-IT qPCR Analysis Kit. To see a list of the available species-specific primers, please visit [www.activemotif.com/chipprimers](http://www.activemotif.com/chipprimers).

#### B. Data Analysis

If the data analysis will not be performed using Active Motif's ChIP-IT qPCR Analysis Kit, two other simplified methods of analysis are provided below. Both methods listed require the generation of a standard curve, containing known amounts of DNA, for each primer pair being used in the experiment.

##### Method 1: Fold enrichment of positive primers relative to negative control primers

1. Produce a standard curve by performing qPCR with your primer set on known DNA quantities of Input DNA (from Section B Step 17) in triplicate. Run three to five samples that are 10-fold dilutions, e.g. 0.005 ng, 0.05 ng, 0.5 ng, 5 ng and 50 ng.
2. Run the ChIP and IgG samples along with the dilution series of the Input DNA standards using both positive control primers (known binding sites) and negative control primers (regions of the genome not bound by your protein of interest).
3. Your qPCR instrument will assign values to each qPCR reaction based on the standard curve. If your machine does not average your triplicate reactions automatically, you will need to calculate these averages.
4. Divide the average value from the positive control primer set by the average value of the negative control primer set to obtain your fold enrichment.

##### Method 2: Express data as a percent of input

1. Produce a standard curve by performing qPCR with your primer set on known DNA quantities of Input DNA (from Section B Step 17) in triplicate. Run three to five samples that are 10-fold dilutions, e.g. 0.005 ng, 0.05 ng, 0.5 ng, 5 ng and 50 ng.

2. Run the ChIP and IgG samples along with the dilution series of the Input DNA standards using both positive control primers (known binding sites) and negative control primers (regions of the genome not bound by your protein of interest).
3. Your qPCR instrument will assign values (in ng) to each qPCR reaction based on the standard curve. If your machine does not average your triplicate reactions automatically, you will need to calculate these averages.
4. For each qPCR reaction you will have used a percentage of your total ChIP DNA. In order to calculate the amount in the whole reaction, divide the elution volume of the entire ChIP reaction by the volume used in the qPCR reaction (*e.g.* if you eluted ChIP DNA in 200  $\mu$ l and used 5  $\mu$ l in the qPCR reaction the formula is  $200/5 = 40$ ). Then, multiply the average qPCR quantity by this number (*e.g.* qPCR quantity in ng  $\times$  40).
5. To express data as a percent of input, divide the adjusted values from Step 4 above by the amount of DNA that went into the ChIP reaction and then multiply by 100%. (*e.g.* if 20  $\mu$ g was used in the ChIP reaction this is equivalent to 20,000 ng of chromatin. The calculation would be the adjusted value from Step 4 divided by 20,000 ng and then multiplied by 100). Typical percent of input recovered values are 0.05% to 1%.



## Section H. Troubleshooting Guide

Problem/question	Recommendation
At what points in the protocol can I stop?	The protocol may be stopped and samples stored at the times and temperatures below: <ol style="list-style-type: none"> <li>1. After formaldehyde fixation and centrifugation (intact cell pellet), -80°C.</li> <li>2. After flash freezing, -80°C.</li> <li>3. After chromatin shearing, -80°C.</li> <li>4. After DNA clean up, -20°C.</li> </ol>
Poor yield of sheared chromatin.	<p>Insufficient cell numbers were used. Repeat chromatin preparation using a larger number of cells.</p> <p>Nuclei not released. Monitor cell lysis under a microscope. Generally, the more cells that are lysed, the higher the sheared chromatin yield. Vortex cells at highest setting for 30 seconds in the presence of Swelling buffer and Detergent.</p> <p>Sonication samples were emulsified. Avoid emulsification by turning up the power of the sonicator gradually. If a chromatin preparation becomes emulsified inadvertently, discontinue shearing and centrifuge the sample for 4 minutes at 8,000 rpm in a 4°C microcentrifuge to remove trapped air.</p> <p>Use fresh formaldehyde when preparing Fixation Solution</p>
Shearing efficiency is not clear from gel analysis.	<p>Material is stuck in the wells, and smears or streaks are seen from the top to bottom of the lane. The sheared chromatin needs to have the cross-links reversed, protein removed (Proteinase K) and RNA removed (RNase), followed by DNA purification.</p> <p>High molecular weight products. Decrease the size of the fragments by re-sonicating the sample. If an alternative reverse cross-linking method was used, or the 20 minute incubation at 100°C in NaCl was omitted prior to running the agarose gel for analysis, please repeat the input chromatin preparation and follow the manual instructions.</p>
Performing ChIP with a large volume of chromatin.	This is not recommended. It is better to set up several small ChIP reactions (240 µl each) and pool the samples during loading of the DNA purification column, rather than trying to ChIP a single large sample. Do not perform a single scaled-up reaction, as the capture efficiency is lower.
ChIP DNA does not turn bright yellow following the addition of 3 M sodium acetate	If the color is light orange or violet, this indicates the pH is too high. Add more 3 M sodium acetate 5 µl at a time, mixing after addition until the color is bright yellow. This step is crucial to the success of DNA binding and purification. For a full color image please see the manual for Active Motif's Chromatin IP DNA Purification Kit Catalog No. 58002 available online at our website <a href="http://www.activemotif.com">www.activemotif.com</a> .
High background.	<p>Chromatin not sheared enough. Shearing should produce DNA fragments that are small enough to exclude background from neighboring chromosomal sequences, but still large enough that there is a good possibility your amplicon remains intact. We recommend 200-1200 bp fragments. If the DNA fragments are too large, the background is increased. Consider increasing the number of pulses for sonication. Check the fragment size on a gel to assess your shearing efficiency.</p> <p>Antibody issue. Too much antibody relative to the amount of chromatin in the ChIP reaction. Excess antibody will result in more non-specific binding, which will be detected as increased background.</p>

Problem/question	Recommendation
<p>Poor or no enrichment with target antibody.</p>	<p>Too little chromatin. Generally, we recommend using 10 - 30 µg of chromatin per ChIP reaction. ChIP reactions should not exceed 50 µg per IP reaction. Be sure to quantitate the concentration of the sheared chromatin sample(s) being ChIP'd to ensure both that adequate chromatin is used per sample, and that equal mass quantities of chromatin are used in each ChIP.</p>
	<p>Antibody is not ChIP validated. The antibody does not efficiently recognize fixed proteins, either because the epitope is destroyed by fixation or because the epitope is masked by other proteins in a larger complex. To assist in ChIP validating an antibody, it is very useful to use a positive control antibody such as Histone H3K4me3 (Catalog No. 39915) and a negative IgG from the same species, and primers that have been proven to work in the type of PCR being used. Active Motif offers species-specific ChIP-IT Control qPCR Kits for antibody validation (Catalog Nos. 53026, 53027 &amp; 53028).</p>
	<p>Low-affinity antibody. Use a different antibody.</p>
	<p>Antibody affinity to protein G is weak. Individual monoclonals have variable binding affinities to protein G, which are pH dependent; the optimal pH may vary for each Ig. For those with low to medium affinity, capture efficiency by protein G can be dramatically improved through use of our Bridging Antibody (Catalog No. 53017). This antibody is a rabbit anti-mouse pAb that recognizes all subclasses of mouse immunoglobulins. If your IgG has a weak/medium affinity to protein A or G, the Bridging Antibody will increase antibody capture by the beads without increasing background.</p>
	<p>Problems with PCR. Check the conditions of the PCR master mix used. Confirm the amplified sequence for the positive control primer set is bound by the antibody target. Identify other binding sites.</p>
<p>No PCR products for the ChIP'd samples (but the Input DNA yields the correct PCR product)</p>	<p>Increase the amount of chromatin used in the ChIP reaction, the amount of antibody used, or both.</p>
	<p>Use a different antibody.</p>
<p>No PCR products with real-time PCR</p>	<p>Confirm the species specificity and efficiency of your primers. You may need to redesign your primers. Primers that work in end point PCR do not always work in qPCR.</p>
	<p>No ethanol in DNA Purification Wash Buffer. Make sure that ethanol has been added to the DNA Purification Wash Buffer prior to first use.</p>

## Section I. Related Products

ChIP-IT® Kits	Format	Catalog No.
ChIP-IT® Express	25 rxns	53008
ChIP-IT® Express Enzymatic	25 rxns	53009
ChIP-IT® Express Shearing Kit	10 rxns	53032
ChIP-IT® Express Enzymatic Shearing Kit	10 rxns	53035
ChIP-IT® High Sensitivity	16 rxns	53040
ChIP-IT® qPCR Analysis Kit	10 rxns	53029
ChIP-IT® ChIP-Seq	10 libraries	53041
ChIP-IT® FFPE	16 rxns	53045
ChIP-IT® FFPE Chromatin Preparation Kit	5 rxns	53030
ChIP-IT® Express HT	96 rxns	53018
Re-ChIP-IT®	25 rxns	53016
RNA ChIP-IT®	25 rxns	53024
Chromatin IP DNA Purification Kit	50 rxns	58002
EpiShear™ Probe Sonicator	110 V	53051
EpiShear™ Cooled Sonication Platform, 1.5 ml	1 platform	53080
ChIP-IT® Protein G Magnetic Beads	25 rxns	53014
Protein G Agarose Columns	30 rxns	53039
Siliconized Tubes, 1.7 ml	25 tubes	53036
ChIP-IT® Control qPCR Kit – Human	5 rxns	53026
ChIP-IT® Control qPCR Kit – Mouse	5 rxns	53027
ChIP-IT® Control qPCR Kit – Rat	5 rxns	53028
ChIP-IT® Control Kit – Human	5 rxns	53010
ChIP-IT® Control Kit – Mouse	5 rxns	53011
ChIP-IT® Control Kit – Rat	5 rxns	53012
Ready-to-ChIP HeLa Chromatin	10 rxns	53015
Ready-to-ChIP Hep G2 Chromatin	10 rxns	53019
Ready-to-ChIP K-562 Chromatin	10 rxns	53020
Ready-to-ChIP NIH/3T3 Chromatin	10 rxns	53021
Bridging Antibody for Mouse IgG	500 µg	53017
Dounce Homogenizer	1 ml	40401
Dounce Homogenizer	15 ml	40415

### ChIP-validated Antibodies

For an up-to-date list of over 125 ChIP-validated antibodies, please visit [www.activemotif.com/chipabs](http://www.activemotif.com/chipabs).

Whole Genome Amplification	Format	Catalog No.
GenoMatrix™ Whole Genome Amplification Kit	1 kit	58001
Co-Immunoprecipitation	Format	Catalog No.
Nuclear Complex Co-IP Kit	50 rxns	54001
Universal Magnetic Co-IP Kit	25 rxns	54002
Modified Histones Array	Format	Catalog No.
MODified™ Histone Peptide Array	1 array	13001
Histone Modification FP Binding Assay	Format	Catalog No.
HiLite™ Histone H3 Methyl-Lys9 / Lys27 FP Binding Assay	1 kit	57001

<b>Histone ELISAs</b>	<b>Format</b>	<b>Catalog No.</b>
Histone H3 monomethyl Lys4 ELISA	1 x 96 rxns	53101
Histone H3 dimethyl Lys4 ELISA	1 x 96 rxns	53112
Histone H3 trimethyl Lys4 ELISA	1 x 96 rxns	53113
Histone H3 acetyl Lys9 ELISA	1 x 96 rxns	53114
Histone H3 dimethyl Lys9 ELISA	1 x 96 rxns	53108
Histone H3 trimethyl Lys9 ELISA	1 x 96 rxns	53109
Histone H3 phospho Ser10 ELISA	1 x 96 rxns	53111
Histone H3 acetyl Lys14 ELISA	1 x 96 rxns	53115
Histone H3 monomethyl Lys27 ELISA	1 x 96 rxns	53104
Histone H3 trimethyl Lys27 ELISA	1 x 96 rxns	53106
Histone H3 phospho Ser28 ELISA	1 x 96 rxns	53100

<b>Histone Purification &amp; Chromatin Assembly</b>	<b>Format</b>	<b>Catalog No.</b>
Histone Purification Kit	10 rxns	40025
Histone Purification Mini Kit	10 rxns	40026
Chromatin Assembly Kit	10 rxns	53500
HeLa Core Histones	36 µg	53501

### **Recombinant Methylated, Acetylated and Phosphorylated Histone Proteins**

For an up-to-date list of Recombinant Histone Proteins, please visit [www.activemotif.com/recombhis](http://www.activemotif.com/recombhis).

<b>Histone Acetyltransferase and Deacetylase Activity</b>	<b>Format</b>	<b>Catalog No.</b>
HAT Assay Kit (Fluorescent)	1 x 96 rxns	56100
Recombinant p300 protein, catalytic domain	5 µg	31205
Recombinant GCN5 protein, active	5 µg	31204
HDAC Assay Kit (Fluorescent)	1 x 96 rxns	56200
HDAC Assay Kit (Colorimetric)	1 x 96 rxns	56210

<b>Histone Demethylase Activity</b>	<b>Format</b>	<b>Catalog No.</b>
Histone Demethylase Assay (Fluorescent)	48 rxns	53200

<b>DNA Methylation</b>	<b>Format</b>	<b>Catalog No.</b>
hMeDIP	10 rxns	55010
MeDIP	10 rxns	55009
MethylDetector™	50 rxns	55001
MethylCollector™	25 rxns	55002
MethylCollector™ Ultra	30 rxns	55005
UnMethylCollector™	30 rxns	55004
Hydroxymethyl Collector™	25 rxns	55013
DNMT Activity / Inhibition Assay	96 rxns	55006
Methylated DNA Standard Kit	3 x 2.5 µg	55008
Fully Methylated Jurkat DNA	10 µg	55003
Jurkat genomic DNA	10 µg	55007

## Technical Services

---

If you need assistance at any time, please call Active Motif Technical Service at one of the numbers listed below.

### Active Motif North America

1914 Palomar Oaks Way, Suite 150  
Carlsbad, CA 92008

USA

Toll Free: 877 222 9543  
Telephone: 760 431 1263  
Fax: 760 431 1351  
E-mail: [tech\\_service@activemotif.com](mailto:tech_service@activemotif.com)

### Active Motif Europe

Avenue Reine Astrid, 92  
B-1310 La Hulpe, Belgium

UK Free Phone: 0800 169 31 47  
France Free Phone: 0800 90 99 79  
Germany Free Phone: 0800 181 99 10  
Telephone: +32 (0)2 653 0001  
Fax: +32 (0)2 653 0050  
E-mail: [eurotech@activemotif.com](mailto:eurotech@activemotif.com)

### Active Motif Japan

Azuma Bldg, 7th Floor  
2-21 Ageba-Cho, Shinjuku-Ku  
Tokyo, 162-0824, Japan

Telephone: +81 3 5225 3638  
Fax: +81 3 5261 8733  
E-mail: [japantech@activemotif.com](mailto:japantech@activemotif.com)

Visit Active Motif on the worldwide web at <http://www.activemotif.com>

At this site:

- Read about who we are, where we are, and what we do
- Review data supporting our products and the latest updates
- Enter your name into our mailing list to receive our catalog, *MotifVations* newsletter and notification of our upcoming products
- Share your ideas and results with us
- View our job opportunities

**Don't forget to bookmark our site for easy reference!**